




Maximum Reliability
in Tablet Production
In-process Control

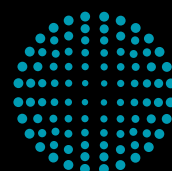

FETTE
COMPACTING
be efficient

The quality requirements for tablet production are constantly increasing. Process equipment from Fette Compacting ensures that your production fully complies with the applicable regulations and effectively reduces the workload of your staff. Extensive automation of all process steps reduces manual intervention and provides for excellent tablet quality.

The Checkmaster makes the automated measurement of the most important tablet parameters possible: weight, thickness, diameter and hardness – and completely integrates them into the in-process control concept.

Your advantages: maximum reliability, the highest quality and efficiency in tablet production.

FEATURES	BENEFIT
+ Fully automatic measurement of the four quality critical parameters: weight, thickness, hardness, diameter	+ Frequent sampling for seamless batch documentation
+ Automatic feedback to tablet weight control loop	+ Easy to clean, rapid availability after a product changeover
+ Precise control of tablet parameters within freely adjustable warning and stop limits	+ Guided validation and calibration process
+ Documented tablet inspection results in your batch report	+ Extremely accurate and reliable weight monitoring
+ 21 CFR Part11 compliant operation via Fette Compacting HMI (Human Machine Interface)	+ All common formats can be tested in standard configuration, customized star-wheels for non-standard formats available
+ GMP-compliant design	



TECHNICAL DATA

Measurement	Weight	Thickness	Diameter	Hardness
Measurement method	Precision weighing cell Sartorius WZ64S	Linear potentiometer	Pressure Load Cell (DMS)	Last Mess Cell (LMZ) with Pressure Load Cell (DMS)
Measuring unit	1 mg / 0.1 mg (adjustable)	mm or inch	mm or inch	N, Kp or Sc
Measuring range	0.01 g – 50 g	1–17 mm	3–18 mm (3–30 mm manuell)	4–400 N standard, optional 4–40 N / 8–800 N
Measuring accuracy	± 1 mg in stable environment, ± 2 mg in unstable environment	± 0.05 mm	± 0.05 mm	0,4–40 N (± 0,1 N) 4–400 N (± 1 N) 8–800N (± 2 N)
Dimensions W x D x H	590 x 400 x 530 mm			
Weight	50 kg (basic)			
Power supply	100–240 V, 50/60 Hz			

4-Parameter Tester, Containment Version

The flexible and modular 4-Parameter Tester is used for fully automated measurement of the four critical parameters: weight, thickness, diameter and hardness.

FEATURES	BENEFIT
+ Separation between the product area and the mechanical and electrical parts	+ CFR21 Part 11 operator-friendly software with reporting, including values / statistics / production tracking
+ Oval, square and rectangular shape centering device	+ New, stable weighing cell – less sensitive to production environments
+ File transfer via Ethernet, USB stick and Bluetooth printing	+ Easy cleaning, validation, calibration and maintenance
+ Local HMI display for parameter modification, fast test review and calibration	+ Extremely quiet test environment – no air turbulence
+ Containment ≤OEB4	



TECHNICAL DATA				
Measurement	Weight	Thickness	Diameter	Hardness
Measurement method	Precision weighing cell Sartorius WZA224-L	Linear Potentiometer	Pressure Load Cell (DMS)	Last Mess Cell (LMZ) with Pressure Load Cell (DMS)
Measuring unit	1 mg / 0.5 mg / 0.1 mg (adjustable)	mm or inch	mm or inch	N
Measuring range	0.01 g – 50 g	3–18 mm	round (3–25 mm) shaped (5–22 mm)	4–400 N or 8–800 N
Measuring accuracy	± 1 mg in stable environment, ± 2 mg in unstable environment	± 0.05 mm	± 0.1 mm	±2 N (800 N ± 4 N)
Dimensions W × D × H	500 × 580 × 990 mm			
Weight	45 kg			
Power supply	100–240 V, 50/60 Hz			

AT50



Tablet testing system that can be fully validated and automatically positions the tablet for measurement of weight, thickness, diameter and hardness.

Integrated Touch-Screen
The AT50 comes standard with an integrated touch screen (EasyTouch™). The EasyTouch™ Software enables easy adjustment of internal testing parameters to adjust to the exact shape of the tablet and speed up the measurement process.

- Main features**
- + Tests can be triggered fully automatically by the tablet press
 - + CFR21 Part11 compliant control via Fette Compacting HMI
 - + IP 52 protective housing
 - + Mettler Toledo weighing cell
 - + Optional rollers for precise tablet positioning on hardness/diameter station
 - + Modular parts exchange
 - + Possible integration into q-doc management software



BENEFIT
+ Automatic measurement of 4 physical parameters
+ Reliable orientation of complexly shaped tablets
+ Full integration with tablet press for self-regulation
+ Easy cleaning and parts exchange
+ Same measurement technology as Sotax laboratory devices
+ Robust design for production environment

TECHNICAL DATA				
Measurement	Weight	Thickness	Diameter	Hardness
Measuring range	0.1 mg – 60 g	2–20 mm	2–25 mm	2–800 N
Resolution	0.1 mg	0.01 mm	0.01 mm	1 N
Dimensions W × D × H	580 × 730 × 527 mm			
Weight	76 kg			
Power supply	100 – 240 V, 50/60 Hz			

NIR Checkmaster

NIR analysis of tablets can significantly cut release times. This technology enables to shorten time and costs on the batch release procedures.

FEATURES

- + Reduction of release times
- + Fully automatic measurement of the critical parameters during the production
- + Tableting is understood and controlled in the sense of the FDA
- + The patented NIRCAL-Software enables a robust and extremely precise calibration

BENEFIT

- + Higher process reliability, test cycles on production process and quality specifications
- + Reduction of production time and costs
- + Shorter release times (Real-Time-Release)
- + Precise positioning for reproducible results



TECHNICAL DATA*

Interferometer type (NIR-Unit)	Polarization interferometer
Detector type	InGaAs, fitted with thermost
Spectral range	12,500–6,000 cm ⁻¹ (recommended 11,520 – 6,000 cm ⁻¹), resp. 800–1,660 nm (recommended 870–1,660 nm)
Resolution	8 cm ⁻¹ (Boxcar-apodization)
Dynamic photometric range	0–6 AU
Wave number accuracy	± 0.2 cm ⁻¹
Typical signal-noise ratio (Rms for spectral segments of 300 cm ⁻¹ in range between 11.000–6.500 cm ⁻¹)	open beam: mean value: 2 × 10 (-5) AU measured with 5 mm spectralon: 10 × 10 (-5) AU
Lamp type	Tungsten-halogen with 2 × 6,000 h working life (double lamp module)
Laser type	HeNe
Dimensions W × D × H	586 × 740 × 581 mm
Weight	107 kg
Power supply	100–240 V, 50/60 Hz

*For the NIR Checkmaster the same technical data as for the Checkmaster apply.

Weightmaster 6.2



Cost-optimized process equipment for in-process control of tablet weight.

FEATURES

- + Fully automatic measurement of tablet weight
- + Automatic feedback to tablet weight control loop
- + Precise control of tablet parameters within freely adjustable warning and stop limits
- + Documented tablet inspection results in your batch report
- + 21 CFR Part11 compliant operation via the Fette Compacting HMI (Human Machine Interface)

BENEFIT

- + Frequent sampling for seamless batch documentation
- + Easy to clean, rapid availability after a product changeover
- + Guided validation and calibration process
- + Extremely accurate and reliable weight monitoring
- + All common formats can be tested in standard configuration, customized star-wheels for non-standard formats available

- + GMP-compliant design

TECHNICAL DATA

Measurement method	Precision weighing cell Sartorius WZ64S
Gauge accuracy	1 mg / 0.5 mg / 0.1 mg (adjustable)
Measuring accuracy	± 1 mg in stable environment ± 2 mg in unstable environment
Measuring range	0.01 g – 50 g
Dimensions W × D × H	480 × 480 × 480 mm
Weight	30 kg
Power supply	100–240 V, 50/60 Hz



Fette Compacting GmbH

Grabauer Strasse 24
21493 Schwarzenbek, Germany
Phone +49 4151 12-0
Fax +49 4151 3797
tablet@fette-compacting.com

Fette Compacting America, Inc.

400 Forge Way
Rockaway N.J. 07866, USA
Phone +1 973 5868722
Fax +1 973 5860450
sales@fetteamerica.com

Fette Compacting America Latina Ltda.

Av. Cambacica, 1200 módulo 10
Parque Imperador
CEP 13097-160
Campinas / SP, Brazil
Phone / Fax +55 19 37969910
contato@fette-compacting.com.br

Fette Compacting Mexico, SA de CV

Adolfo Prieto No. 1638
Colonia Del Valle Sur
03100 Mexico, DF, Mexico
Phone +52 55 40000653
tablet@fette-compacting.com

Fette Compacting (China) Co., Ltd.

No. 9 Shengtong Road,
Moling Sub-District,
Jiangning Development Zone,
211111 Nanjing
Jiangsu Province, P.R.C., China
Phone +86 25 52121818
Fax +86 25 52129951
fcn@fette-compacting.com

Fette Compacting Machinery India Private Limited

A - 406 /407, 4th floor, Atrium 215,
Next to Hotel Courtyard Marriott,
Near J. B. Nagar Metro station,
Andheri – Kurla Road, J. B. Nagar,
Andheri (East)
400 093 Mumbai, India
Phone +91 83 26750355
sales.in@fette-compacting.com

Competence Centre
Plot No S 115, Phase III B
Verna Industrial Estate
Verna, Goa 403 722, India
Phone +91 83 26750355

Fette Compacting Asia Pacific Pte Ltd.

107 Eunos Avenue 3, #01-01
Singapore 409837, Singapore
Phone +65 659 25654
Fax +65 654 71939
infoasiapacific@fette-compacting.com

Fette Compacting Ibérica SL

Avenida Labradores, 1
2ª Planta, Oficina 3
28760 Tres Cantos, Spain
Phone +34 91 8039689
Fax +34 91 3483052
fcib@fette-compacting.com

Fette Compacting France

1, Rue du Centre
93160 Noisy Le Grand, France
Phone +33 155 812121
Fax +33 155 812120
fcf@fette-compacting.com

Fette Compacting Belgium BVBA

Schaliënhoevedreef 1b
2800 Mechelen, Belgium
Phone +32 15 684260
Fax +32 15 684269
fcbe@fette-compacting.com

EuroPharma Machinery Ltd

Unit 12 Highview
Bordon, Hampshire, GU35 0AX
United Kingdom
Phone +44 1420 473344
Fax +44 1420 488030
admin@europharma.co.uk


Fette Compacting Middle East FZE

Jebel Ali Free Zone, Jafza
Lobby 14, Office 308, Dubai
United Arab Emirates
Phone +971 4 8808226
dubai@fette-compacting.com

Fette Compacting North-West Africa

Tour d'Affaire Nord, Centre Commercial
Bab Ezzouar – Regus 4ème Etg.
16000 Algiers, Algeria
Phone +213 770 117 007
fcnwa@fette-compacting.com

www.fette-compacting.com

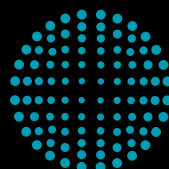

**FETTE
COMPACTING**
be efficient